Isotemp Hybridization Incubators

Small: 13-247-10 (120V), 13-247-11 (100V),
13-247-12 (230V CE), 13-247-13 (230V UK)

Medium: 13-247-20 (120V), 13-247-21 (100V),
13-247-22 (230V CE), 13-247-23 (230V UK)

Dual: 13-247-30 (120V), 13-247-31 (100V),
13-247-32 (230V CE), 13-247-33 (230V UK)

Table of Contents

Information	3
Contact Information	3
Overview	3
Equipment Summary	3
Models Available	4
Rotisseries Available	4
General Specifications	5
Safety Information	6
Unpacking and Installation	9
Unpacking	9
Packing List	9
Location	9
Installation Procedure	10
Operation	11
Front Panel Controls and Indicators	11
Rear Panel Nomenclature	12
General Operation Procedure	13
Power Up	14
Power Up Messages	14
Setting the Temperature	15
Setting the Keypad Protection	15
Setting the Offset Temperature	16
Setting and Reading Temperature Deviation	16
Setting Temperature Deviation On or Off	16
Reading and Resetting Temperature Deviation	17
Preparing the Rotisserie	18
Using the Rotisserie	18
Installing the Rotisserie into the Incubator	19
Loading the Rotisserie	19
Setting the Rotisserie for Variable Axis	20
Unloading the Rotisserie	20
Using the Rocker	21
Installing the Rocker	21
Loading the Rocker	22
Unloading the Rocker	22
Maintenance	23
Cleaning	23
Service	24
Replacement Parts	25
Replacing the Fuses	26
Tightening Loose Belt Tension	26
Fisher Chemical BioReagents	28
Accessories	29

Introduction

This manual provides information and procedures on how to unpack, set up, and operate the Fisher Isotemp brand of Hybridization Incubators.

This manual is intended to be used by Fisher customers responsible for operating the hybridization incubator.

Contact Information

Fisher Scientific, 2555 Kerper Blvd. Dubuque, IA 52001

Technical assistance: 1-800-943-2006
(for guidance in using the product)

Instrument service: 1-800-395-5442
(for repairs and replacement parts)

Customer service: 1-800-766-7000
(to order products and accessories)

Copyright 2000, Fisher Scientific. Fisher reserves the right to change specifications or designs at any time without incurring obligation.

Overview

Equipment Summary

Fisher Isotemp Hybridization Incubators are designed to process nucleic and protein blots derived from hybridization procedures. Incubator heat is accurately controlled to precisely maintain the chamber temperature. Temperature can be set between slightly above ambient and 85°C and is easily set by simple controls on the front panel. A large, bright display shows actual temperature during operation. The incubator features keypad protection, which prevents the set temperature from being changed. This is useful when setting up a processing program where the unit will be unattended. An offset temperature can be programmed to compensate for the actual temperature inside a typical sample loaded into the tubes. The incubator can monitor and store the maximum deviation of the actual temperature from the set temperature. Temperature deviation can be monitored and saved during processing and recalled

INTRODUCTION

later. A separate safety thermostat limits the incubator temperature to approximately 95°C in the event of a temperature controller failure. An indicator on the front panel shows when this condition exists. A rotisserie holds the sample tubes and rotates to provide membrane coverage. Rotisserie rotation speed can be varied by a knob on the front panel. A jog control permits rotating the rotisserie in steps to insert and remove tubes. The incubator has a glass door for observing samples during processing. Interior surfaces are stainless steel for high durability and low maintenance. Small and medium same size units are stackable with an optional stacking plate. The small incubator can be mounted under a cabinet or counter with the optional mounting bracket.

Models Available

Size	Incubator Catalog Number				Sample Tube Capacity		
	at line voltage				with supplied rotisserie		with optional rotisserie
	120 VAC	100 VAC	230 VAC Cont. European	230 VAC UK	35x300 mm	35x150 mm	70x300 mm
Small	13-247-10	13-247-11	13-247-12	13-247-13	4	8	2
Medium	13-247-20	13-247-21	13-247-22	13-247-23	10	20	4
Dual	13-247-30	13-247-31	13-247-32	13-247-33	20	40	8

Rotisseries Available

Different size rotisseries accept different sample tube capacities:

- With the small unit, the supplied rotisserie accepts 4 of the 35x300 mm sample tubes or 8 of the 35x150 mm tubes. An optional rotisserie accepts 2 of the 70x300 mm tubes.
- With the medium unit, the supplied rotisserie accepts 10 of the 35x300 mm tubes or 20 of the 35x150 mm tubes per incubator. An optional rotisserie accepts 4 of the 70x300 mm tubes per incubator.
- The dual unit, accepts two rotisseries, identical to those describe for use in the medium unit.

As an alternative to rotisseries, a rocker is available for agitating samples. The rocker is an option on the medium and dual units and it replaces the rotisserie.

General Specifications

Specifications

Power Requirements: 100 VAC \pm 10%, 50/60 Hz, 120 VAC \pm 10%, 60 Hz, 230 VAC \pm 10%, 50/60 Hz

Small/medium unit: 3.4 A @ 120 VAC, 1.8 A @ 230 VAC

Dual unit: 6.8 A @ 120 VAC, 3.6 A @ 230 VAC

Ambient conditions: +16 to 40°C, up to 90% relative humidity, non-condensing

Temperature range: Ambient +7 to 85°C

Average uniformity: \pm 0.5°C at 55°C within the chamber; \pm 0.25°C at 55°C sample to sample

Heat-up time: 20 minutes from ambient to 65°C

Rotisserie speed: 5-15 rpm

Maximum altitude: 2000 m

Over-voltage category: II (IEC 664)

Pollution degree: 2 (IEC 664)

Exterior dimensions:

W x L x H (nominal) Small: 18 x 16.5 x 17 inches (46 x 42 x 44 cm)
Medium: 18 x 20 x 20.5 inches (46 x 51 x 53 cm)
Dual: 18 x 20 x 41 inches (46 x 51 x 104 cm)

Weight: Small: 45 lbs (21 kg)
Medium: 55 lbs (25 kg)
Dual: 105 lbs (48 kg)

Note: (230 VAC units only)

CE products meet the relevant EC harmonized standards for safety (IEC1010-1/EN61010) and EMC (EN61326-1).

Safety Information

The following symbols may be used on the device:

Indicates that user must read and understand the instructions.

Indicates that surfaces may be hot.

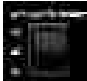
Indicates a shock hazard from dangerous voltage.

Indicates a pinch hazard may exist.

The following labels are used on the device:

	<u>Label</u>	<u>Appearance</u>	<u>Function</u>
Controls	Power		Controls line power to the unit.
	Rotisserie Motor		Controls the motion of the rotisserie motor.
	Rotisserie Speed		Controls the rotisserie's speed of rotation.
	Temperature Control		These controls are associated with the heating function. Used with temperature display (see below).
Indicators	Temperature Display		Indicates the temperature of the incubator chamber. Used with temperature controls (see above). NOTE: The red bar to the left of the display indicates a negative value. Used in calibrating the unit.
	Alarm	Alarm	Indicates that the unit is operating in the alarm condition (the indication is only seen during the alarm condition).

These instructions contain important operating and safety information. The user must carefully read and understand these instructions before using the hybridization incubator.

The hybridization incubator has been designed to optimize function, reliability, safety, and ease of use. It is the user's responsibility to install and use the hybridization incubator properly. Follow these safety instructions regarding the hybridization incubator:

- The hybridization incubator must be used as described in this manual. If used in a manner other than described in this manual, protection provided by the device may be impaired.
- User must completely read, understand, and follow the instructions in this manual before operating the hybridization incubator.
- This equipment is for indoor use only.
- If you plan to use the unit with any procedure that will give off fumes, be sure to operate the unit in a fume hood or with proper ventilation.
- Connect the unit to a properly grounded electrical outlet of the correct voltage and current handling capacity. Check the nameplate on the back of the unit for the voltage and current rating.
- Do not remove or modify the grounded power plug. Use only properly grounded outlets to avoid a shock hazard.
- Disconnect the unit from the power outlet before performing any cleaning, maintenance, or service.
- The hybridization incubator is not intended for use with flammable materials. If used with flammable materials, fire or explosion hazard may result.
- If radioactive materials are used in processing, user must comply with all regulations regarding the acquisition, use, and disposal of such materials.

SAFETY INFORMATION

- Do not continue to use the incubator if the temperature controller fails, as evidenced by the red ALARM indicator lighting.
- Use appropriate hand and eye protection when handling hazardous chemicals.
- Rotisserie and shaker accessory may pose potential pinch hazard. Always turn rotisserie motor off prior to handling rotisserie or shaker.
- The interior of the incubator and its contents can reach temperatures that can cause burns. Avoid contact. The unit can remain hot without visual indication for some time after power is switched off.

CAUTION! DOUBLE POLE NEUTRAL FUSING.

- The power entry module located on the rear of the unit contains two fuses, one located in each side of the AC line.
- To ensure stability of Dual Models (13-247-30, -31, -32 and -33), secure unit to a wall using the included wall bracket.
- Use caution when moving incubator. Note weight of unit to be moved and obtain help if necessary. Tilt incubator to allow grasping under bottom of unit at the sides or back. Do not lift by grasping under plastic bezel at front of unit. Dual unit may require a third person to prevent unit from tipping while being moved.

Unpacking and Installation

Unpacking

Each hybridization incubator is shipped in a single carton. When unpacking the unit, check each item against the packing list below. If something is missing, contact Fisher customer service at 1-800-766-7000, identifying the part by name and catalog number.

Part numbers are listed in the Replacement Parts and Accessory Items sections located at the end of this manual.

Packing List

Small	Quantity		Item supplied
	Medium	Dual	
1	1	1	Hybridization incubator
1	1	2	Drip tray
1	1	2	Rotisserie
1	1	1	35mm x 150mm hybridization tube (Cat. No. 13-247-150)
1	1	1	35mm x 300mm hybridization tube (Cat. No. 13-247-300)
2	2	4	Spare tube clips
1	1	1	Power cord
1	1	1	Instructions
1	1	1	Warranty card
0	0	1	Wall Bracket

Location


Follow these guidelines when choosing a location for the hybridization incubator. The location must:

- Provide an adequate source of power. Check the label on the back of the unit for voltage and current requirements.
- Provide minimum air clearance as follows: 0.5 inch (13 mm) from top, 2 inches (51 mm) from back, and 2 inches (51 mm) on each side.
- Provide adequate clearance for loading and unloading samples.


UNPACKING AND INSTALLATION

- Be level and capable of supporting the weight of the unit when loaded with samples.
- Be free of drafts and wide ambient temperature variations such as near a heater or air conditioning vents.
- Be near sufficient counter space for sample preparation, loading, and unloading.
- Provide a fume hood if hazardous fumes are anticipated when using the hybridization incubator.

Figure 1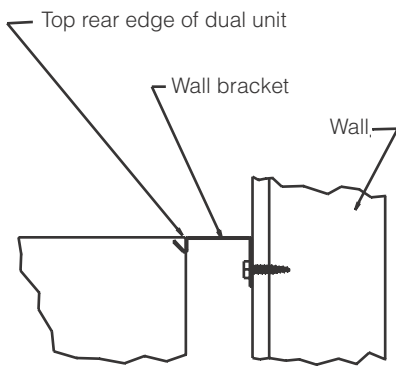

OR

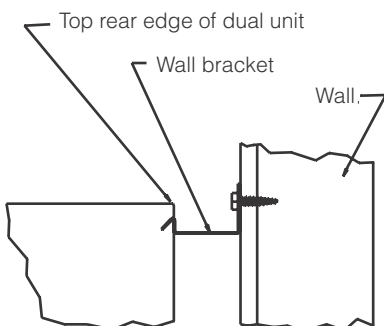
Installation Procedure

1. Carefully unpack the unit and check contents against the packing list.
2. Remove all packaging material before using the incubator.
3. Place incubator on a stable, level surface. To ensure stability of Dual Models (13-247-30, -31, -32, -33), secure unit to a wall using the included wall bracket.
4. Install rotisserie or rocker.
5. Install drip tray inside incubator chamber.
6. Plug the power cord into a suitable grounded electrical outlet.
7. Check out system operation before using unit.
8. Complete and return the warranty card.

Operation

Front Panel Controls and Indicators

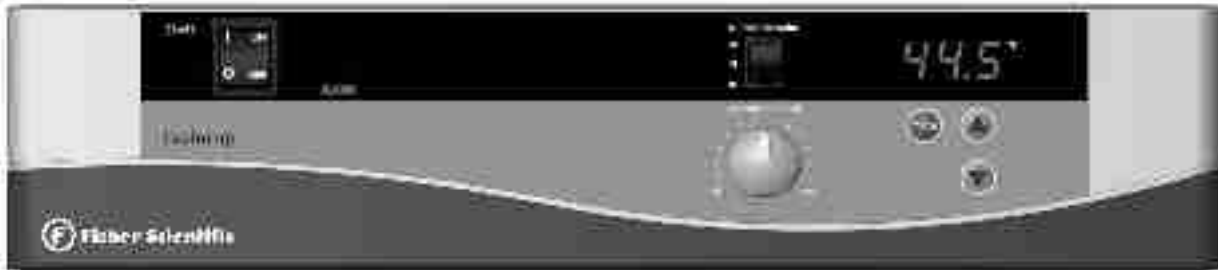Control or indicator

What it does

How to use

power switch

Controls line power to unit.

1 = power on
0 = power off

rotisserie motor switch

Controls motion of rotisserie

off = stops rotisserie
on = runs rotisserie at the speed set on rotisserie speed knob
jog = moves rotisserie a fraction of a turn each time switch is pressed.

rotisserie speed knob

Controls speed of rotisserie

Turn knob to the desired speed for the rotisserie. For slower speeds, turn the knob counterclockwise toward **min.** For faster speeds, turn the knob clockwise toward **max.**

temperature display

Shows actual or set temperature, depending on actions taken on other keys.

During power up, displays actual temperature of incubator. If you press either of the arrow keys, shows set temperature.

Shows messages to indicate menu functions.

Use in conjunction with menu key to access other functions.

Shows status and error messages. Messages)

Take action depending on message. (See Power Up

Front Panel Controls and Indicators (cont.)

Control or indicator
temperature control
menu key

What it does
Advances the display through
through the functions.

How to use
Press the menu key to step all functions.
While each message is displayed, you can
set a value with the arrow keys.

CAL = Temperature offset

The entered value is added to the
measured temperature to allow calibrating
the sample temperature.

Pro = Keypad protection

Choose on or off. When on, set
temperature cannot be changed. When
off, set temperature can be changed.

dEn = Deviation enable

When on, the system stores the maximum
temperature deviation from the set
temperature. When off, the deviation will
not be stored.

dE = Maximum deviation from set
temperature since the last time the
value was reset.

When deviation is enabled (see above),
display shows the maximum deviation
temperature since this function was last
enabled. If deviation enable is set to off,
this function is skipped. Pressing either
arrow key while viewing the deviation
resets the value to zero.

temperature control
arrow keys

Display or change the set temperature

Press either arrow key once to display the
set temperature for two seconds.

Hold the up arrow key to increase the set
temperature. Hold the down arrow key to
decrease the set temperature.

ALARM indicator

Alarm

Shows when there is an alarm
condition

If the ALARM indicator lights, switch off
power and call Fisher Service. Primary
heater control has failed and unit is
operating on safety thermostat.
Temperature is approximately 95°C.

Rear Panel Nomenclature

Item
power entry module/
line filter/fuse holder

What it does
Accepts power cord for providing
line power to unit. Filters incoming
and outgoing power line noise. Holds
fuses for over-current protection.

How to use
Plug mating end of power cord into this
connector. Plug other end into suitable
power outlet.

General Operation Procedure

Follow this general procedure when operating the hybridization incubator:

1. Read, understand, and follow all safety information.
2. Install rotisserie or rocker. See “Installing the Rotisserie into the Incubator” and “Rocker Installation” in the Operation section.
3. Install drip tray and leave in place at all times.
4. Power up system. See “Power Up” in the Operation section.
5. Set temperature as desired for processing and allow incubator to reach and stabilize at desired temperature. See “Setting the Temperature” and “Setting the Temperature Offset” in the Operation section.
6. Load samples into rotisserie or rocker. See “Loading the Rotisserie” in the Operation section. To set the rotisserie for variable axis, see details under “Using the Rotisserie” and “Loading the Rocker.”
7. Run processing procedure as desired. If you want to set keypad protection, see “Setting the Keypad Protection” under the Operation section. For details on temperature deviation, see “Setting and Reading Temperature Deviation” in the Operation section.
8. At conclusion of processing, unload samples from rotisserie or rocker. For details on rotisserie and rocker unloading, see “Unloading the Rotisserie” and “Unloading the Rocker” in the Operation section.
9. Clean incubator for next use. See “Cleaning the Unit” in the Operation section.

Power Up

Follow this procedure to power-up the incubator.

1. Set the power switch to the I position. Display alternates between **oFF** and the measured temperature. If you get any other message, see “Power Up Messages” below.
2. Press any key to stop the **oFF** display. The system displays the actual temperature inside the incubator.

Power Up Messages

When you switch on the power to the hybridization incubator, the display shows one of these messages. Also, if the power ever goes off while the incubator is operating, one of these messages will be displayed:

Message	What it means	What you should do
Display alternates between oFF and the measured temperature.	Power was applied after system had been switched off. If message appears during operation, indicates that power went off.	Press any key to stop the oFF display. Incubator continues operating normally. This is the normal power-up message.
Display alternates between oFF and Alr (alarm).	Measured temperature is more than 3.0° C above the set temperature. May appear if set temperature is changed to a value that is more than 3.0° C below the actual temperature.	Press any key to stop the oFF display. Display alternates between Alr and the actual temperature. Allow incubator to cool. If problem is not the result of setting the temperature lower, have incubator serviced.
Display alternates between oFF and HI .	Measured temperature is more than 99.9° C.	Press any key to stop the oFF display. System will not operate and will continue to display HI . Allow incubator to cool. If problem continues, have incubator serviced.
Display alternates between oFF and tc (temperature control).	There is a problem with either the temperature sensor or the incubator controller.	Press any key to stop the oFF display. System will not operate and will continue to display tc . If problem continues, have incubator serviced.

Setting the Temperature

1. Press and release either of the arrow keys to display the set temperature for two seconds.
2. With the set temperature displayed, to increase the set temperature, press and hold the up arrow for at least two seconds. To decrease the set temperature, press the down arrow. Temperature can be set between 0.0 and 85.0°C.
3. Release the arrow keys when the temperature is set as desired. After ten seconds, the display shows the actual temperature. To return to the actual temperature immediately, press the menu key.

Setting the Keypad Protection

When keypad protection is on, the arrow keys are disabled and the set temperature cannot be changed. This is useful when setting up a processing program where the unit will be unattended. Follow this procedure to set keypad protection on or off:

1. Press the menu key twice so the display shows **Pro** (protection).
 - If keypad protection is on, display shows **on**.
 - If keypad protection is off, display shows **oFF**.
2. To set keypad protection on or off, press either the up or down arrow key.
 - Display shows the new status of keypad protection.
 - After ten seconds, the display returns to the actual temperature.
 - To return to the actual temperature immediately, press the **menu** key.

Setting Offset Temperature

The incubator has been calibrated to display the actual temperature at the center of the chamber. It has been calibrated at 55.0°C therefore the display will be most accurate at this temperature setting. You may wish to recalibrate the incubator to obtain the highest accuracy at another temperature, or to set the temperature to read the actual temperature inside a typical sample inside a sample tube. To do this, use a thermometer to determine the temperature in the chamber or inside the sample, then follow this procedure to set the offset.

1. Note the difference between the displayed temperature and the sample temperature. For example, if the displayed temperature is 45.0°C and the sample temperature is 42.0°C, the offset is -3.0°C. This means that the display should read 3.0°C lower than it now shows.
2. Press the menu key once so the display shows **CAL**.
3. Press and hold either the up or down arrow key to set the desired offset. In the example above, you would set the offset to -3.0°C.
4. Check the temperature accuracy again and change the offset as necessary to achieve the desired temperature.

Setting and Reading Temperature Deviation

With the temperature deviation mode on, the maximum deviation of the actual temperature from the set temperature will be stored and displayed when recalled. With this function off, temperature deviation will not be displayed.

Setting Temperature Deviation On or Off

Follow this procedure to set temperature deviation on or off before processing. After setting on, the maximum tem-

perature deviation will be stored (even with power switched off) until this function is set off or the value is reset.

1. Press the menu key three times so the display shows **dEn** (deviation enable).
 - If temperature deviation is on, display shows **on**.
 - If temperature deviation is off, display shows **oFF**.
2. To set temperature deviation on or off, press either the up or down arrow key.
 - Display shows the new status of temperature deviation.
 - After ten seconds, the display returns to the actual temperature.
 - To return to the actual temperature immediately, press the **menu** key.

Reading and Resetting Temperature Deviation

If temperature deviation is set on, follow this procedure to read the maximum temperature deviation since this function was enabled. You can also follow this procedure to reset the function to start monitoring the temperature deviation again.

1. Press the menu key four times so the display shows **dE** (temperature deviation).

Display shows the maximum temperature deviation from the set temperature since the last time the value was reset.

2. To reset the temperature deviation monitoring function, press either the up or down arrow key while viewing the deviation value.

Display shows temperature deviation of **0.0**.

OPERATION

After ten seconds, the display returns to the actual temperature.

To return to the actual temperature immediately, press the **menu** key.

Using the Rotisserie

Preparing the Rotisserie

To install the tube holders to the rotor discs:

1. Lift tab and slide tube holder onto rotor disc until the tube holder clicks into place. See Figures 2 and 3 below.

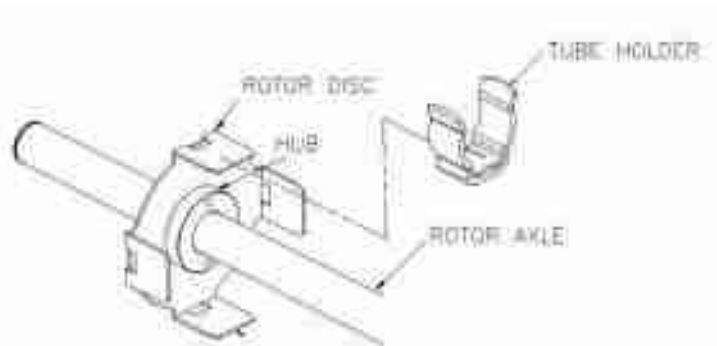
Figure 2

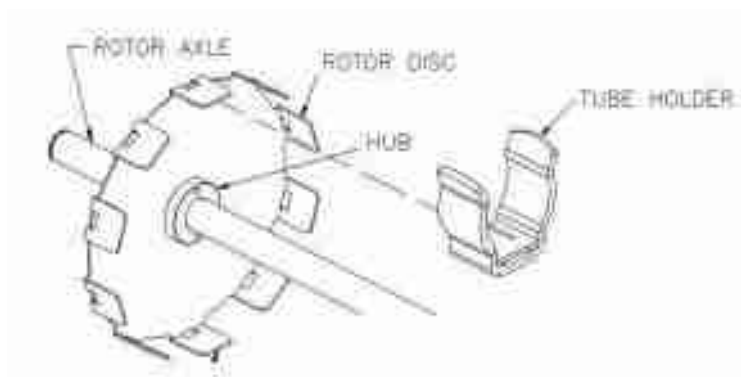
Figure 3

Figure 4

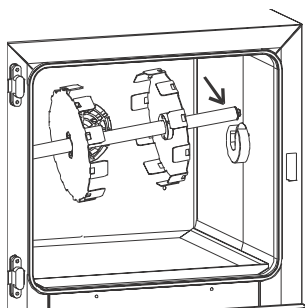
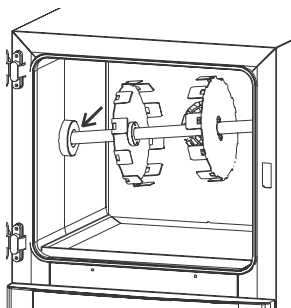
Figure 5

Installing the Rotisserie into the Incubator

Follow this procedure to install a carousel into the incubator.

1. Set the power switch off and open the incubator door.
2. Place the rotisserie into the incubator with the smooth shaft end on the left, gear on the right.
3. Line up the left side of the shaft with the pivot point, then push the rotisserie to the left against the spring, as shown in Figure 4.
4. Line up the right side of the shaft with the gear drive socket, as shown in Figure 5. Release the rotisserie and allow the shaft to be pushed into the socket.

Loading the Rotisserie

With the empty rotisserie in the incubator load the sample tubes as follows:

1. Set the power switch on, set the rotisserie motor switch off, and open the incubator door.
2. Place a sample tube against the rotisserie clips facing the front of the unit. If you need to position the rotisserie, press the rotisserie motor switch to the jog position.
3. Center the sample tube from left to right, then press the tube into the clips until the tube snaps in place.
4. Use the rotisserie motor switch to jog the rotisserie so that an empty position faces the front of the unit.
5. Install another sample tube the same way. Center the tube from left to right, then press the tube into the clips until the tube snaps in place.
6. Repeat jogging the rotisserie and inserting tubes until all of the sample tubes are installed.

Note


Use the rotisserie speed knob to change the jog speed.

Don't manually rotate the rotisserie. Use the jog position of the rotisserie motor switch.

For best operation, tubes should be loaded symmetrically to balance the rotisserie.

OPERATION

7. Close the incubator door, set the temperature, and set the rotisserie motor switch on to start processing.
8. At the conclusion of processing, set the power switch off, open the incubator door, and unload the rotisserie.

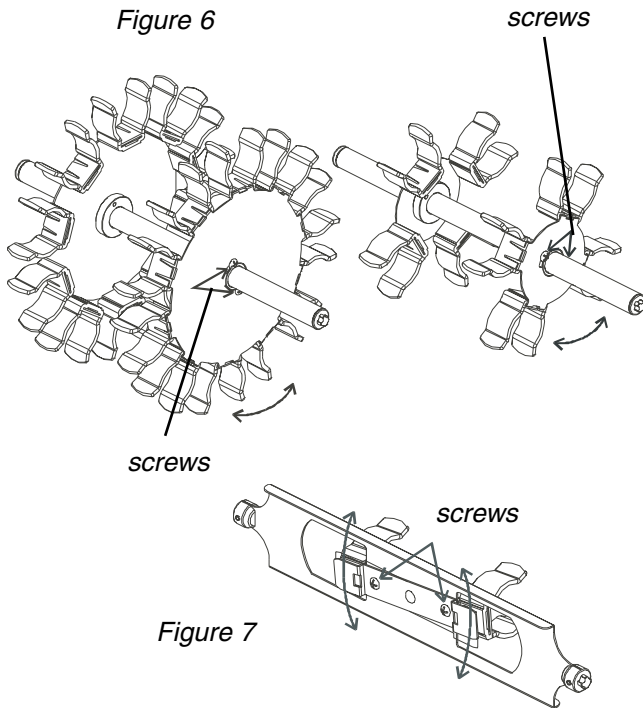Setting the Rotisserie for Variable Axis

With lower volumes in the sample tubes, you may wish to set the rotisserie for variable axis motion to assist in hybridization. Follow this procedure to change the rotisserie motion from straight motion, where the sample tubes remain level, to variable axis motion, where the sample tubes are at an angle.

1. Set the power switch off, then remove the rotisserie from the incubator.
2. Using a Phillips screwdriver, loosen the two screws on one or both ends of the rotisserie hubs, depending on the desired offset. These screws are shown at left.

If you have the optional rotisserie for the small incubator, loosen the two screws near the middle of the rotisserie, shown at left.

3. Set the rotisserie to the desired offset. You may want to install a sample tube and move the rotisserie to check the setting.
4. Tighten the screws so the rotisserie does not twist.
5. Install the rotisserie into the incubator.

Unloading the Rotisserie

After running the program, follow this procedure to unload the rotisserie:

1. Open the incubator door. Have power switched on, and the rotisserie motor switched off.

Note

Don't manually rotate the rotisserie. Use the jog switch.

2. Use the rotisserie motor switch to jog the rotisserie so that a sample tube faces the front of the unit.
3. Pull the sample tube straight out of the rotisserie.
4. Again use the rotisserie motor switch to jog the rotisserie so that a sample tube faces the front of the unit then remove the tube from the rotisserie.
5. Repeat jogging the rotisserie and removing tubes until all of the sample tubes are removed.
6. Set the power switch off.
7. Clean the unit in preparation for the next use.

Using the Rocker

As an alternative to rotisseries, a rocker is available for agitating samples. The rocker is an option on the medium and dual units and it replaces the normal rotisserie.

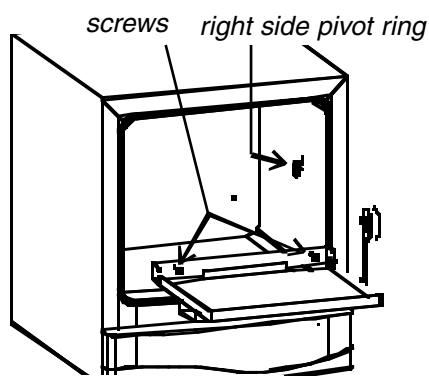
Figure 8

Installing the Rocker

Follow this procedure to install the rocker:

1. Set power switch off and open the incubator door.
2. Remove the rotisserie from the incubator: Push the rotisserie to the left, lift the right side out of the socket, then move the rotisserie slightly to the right and lift it out of the incubator.
3. Place the rocker in the incubator and secure it to the back of the incubator with the two existing screws from the back panel, as shown in Figure 8.
4. Remove the right side pivot ring from inside the incubator: Turn the ring a fraction of a turn counterclockwise, then pull the ring away from the side of the incubator.

OPERATION

Note

When using the rocker, the total sample weight should not exceed 7 lbs (3.2 kg). Distribute the weight evenly, balanced, and centered.

5. Install the rocker pivot ring: Line up the pivot ring spline with the exposed socket then press in and tighten the center fastener.
6. Set power switch on and set the rotisserie motor switch on. Make sure the rocker functions properly with no interference.

Loading the Rocker

With the empty rocker in the incubator, load the samples as follows:

1. Set the power switch on, set the rotisserie motor switch off, and open the incubator door.
2. Install samples into the rocker. If you need to move the rocker for easier loading, use the rotisserie motor switch to jog the rocker.
3. Close the incubator door, set the temperature, and set the rotisserie motor switch on to start processing.

Unloading the Rocker

Follow this procedure to unload the rocker:

1. Open the incubator door. Have power switched on, and the rotisserie motor switched off.
2. Use the rotisserie motor switch to jog the rocker if necessary to assist in removing samples.
3. Set the power switch off.
4. Clean the unit in preparation for the next use.

Maintenance

Cleaning

All of the incubator chamber walls are stainless steel. Clean the interior of the incubator regularly to keep the stainless steel from corroding.

The exterior of the unit is painted metal and plastic. Clean the exterior with warm water or a mild cleaning solution. Follow these guidelines when cleaning the unit:

- Unplug the unit, leave the door open, and allow the incubator to cool before cleaning.
- Remove the drip tray from inside the chamber.
- Clean all surfaces and the drip tray with mild detergent (such as dishwashing detergent) and warm water.
- If radioactive materials have been used, clean the incubator using an appropriate cleaner in an approved manner.
- Do not scrub any surface with steel wool. Steel wool leaves small metal particles behind that will rust, causing the surface to look rusty.
- Where there is a heavy coating of residue on the stainless surfaces, or where there is evidence that corrosion is beginning, use a stainless steel cleaner.
- Rinse and wipe all surfaces with clean water and allow to air dry.

Service

If you need replacement parts or repairs on your hybridization incubator, call the Fisher Service Division at 1-800-395-5442. We suggest that you refer any servicing needs for your hybridization incubator to qualified personnel.

If you are having problems using the incubator, follow these general procedures to track down the cause.

Symptom	Possible Causes of Problem
No power	Unit not plugged in; fuse(s) blown; no power at the outlet; incorrect power (make sure outlet matches label on back of unit); defective power cord; defective power switch.
No heating or heating up slower than normal	Defective heater coil; defective temperature controller board.
Fan doesn't run	Defective fan motor.
Rotisserie doesn't turn	Rotisserie motor switch set to off; defective rotisserie motor; broken drive belt; defective motor controller board.
Display alternates between actual temperature and Alr (alarm).	Temperature setting has been reduced below current chamber temperature. Defective solid-state relay.
ALARM light on	Defective thermocouple; defective temperature controller board; defective solid state relay, defective safety thermostat.

Replacement Parts

Use only genuine Fisher parts in your hybridization incubator. When ordering parts from Fisher, use these service part numbers. To order parts, call Fisher at 1-800-766-7000.

Item	Service Part No.
Temperature controller board (100 V and 120 V units)	SPN103703
Temperature controller board (230 V units)	SPN104627
Motor controller board	SPN103706
Solid state relay	SPN88616
Heater, 100/120 VAC	SPN103644
Heater, 230 VAC	SPN105019
Rotisserie motor	SPN103686
Timing belt	SPN103977
Fan motor, 100/120 VAC	SPN103990
Fan motor, 230 VAC	SPN103991
Temperature sensor	SPN95603
Safety thermostat	SPN103986
Power switch	SPN102629
ALARM light, 100/120 VAC	SPN30244
ALARM light, 230 VAC	SPN23358
Rotisserie motor switch	SPN103989
Rotisserie speed potentiometer	SPN103987
Knob for rotisserie speed control	SPN103867
Power entry module, (for all 120V units except on 13-247-30 and -31)	SPN102486
Power entry module, (for 13-247-30 and -31)	SPN104055
Line Filter (for 13-247-12, -13, -22, -23, -32, and -33)	SPN104579
Fuse, 5A, 250V, 5 x 20mm, fast acting	SPN102487
Circuit Breaker, 8A, (for 13-247-30 and -31)	SPN103988
Power cord, North American	SPN40521
Power cord, Euro	SPN83115
Power cord, UK	SPN83117
Door gasket	SPN103967
Front panel overlay, with "wave"	SPN103974
Front panel overlay, no "wave"	SPN103975
Rotisserie tube clip, 35 mm	SPN103958
Rotisserie tube clip, 70 mm	SPN104008

Warning

Double pole neutral fusing.

Replacing the Fuses

There are two fuses located in the power entry mode on the rear of most units. To replace fuses:

1. Turn the power switch to the **OFF** position and pull the power cord out of the module.
2. Fit a thin blade screwdriver under the bottom portion of the fuse holder.
3. Using a screwdriver, slide the fuse holder out of the module.
4. Replace the bad fuse(s).
5. Slide the fuse holder back into the module until it snaps into place.

Tightening Loose Belt Tension

For Cat # 13-247-10, 11, 12 and 13, perform the following:

1. Turn the power switch to the OFF position and pull the power cord out of the module.
2. Lay the incubator on its right side.
3. Using a screwdriver, remove the screws, then remove the chassis bottom.
4. Using a 11/32 nut driver, loosen the two nuts that hold motor bracket to the chassis.
5. Slide the motor to the rear to tighten belt tension, then tighten nuts.
6. Replace chassis bottom.

For Cat # 13-247-20, 21, 22, 23, 30, 31, 32 and 33, perform the following;

1. Turn the power switch to the OFF position and pull the power cord out of the module.
2. View the incubator from the rear.

3. Using a screwdriver, remove the screws, then remove the chassis back cover.
4. Using a 11/32 nut driver, loosen the two nuts that hold motor bracket to the chassis.
5. Slide the motor to the rear to tighten belt tension, then tighten nuts.
6. Replace chassis back cover.

Fisher Chemical BioReagents

The following BioReagents may be used with nucleic acid hybridizations. These solutions are all guaranteed RNase, DNase, and Protease free. They are sterilized through 0.2 um filter and/or autoclaved, and they are manufactured under ISO 9002 guidelines. When ordering these materials from Fisher, use the catalog numbers below. To order these materials, call Fisher at 1-800-766-7000.

Item	Size	Catalog No.
Denhardt's Reagent	150 mg	BP520-5
Denhardt's Reagent, 50X	5 ml	BP515-5
Dextran Sulfate, Sodium Salt	100 g	BP1585-100
	500 g	BP1585-500
Formamide	25 ml	BP531-25
	500 ml	BP531-500
Hybridization Cocktail, 50% Formamide	50 ml	BP1575-50
Polyvinylpyrrolidone	100 g	BP431-100
Bovine Serum Albumin (Fraction V)	100 g	BP1605-100
Sodium Dodecyl Sulfate, 20% Solution (SDS)	200 ml	BP1311-200
	1 liter	BP1311-1
Saline Sodium Citrate, 20X Solution (SSC)	1 liter	BP1325-1
	20 liter	BP1325-20
Saline Sodium Phosphate-EDTA, 20X Solution (SSPE)	1 liter	BP1328-1
	20 liter	BP1328-20
Tris Borate EDTA, 10X Solution (TBE)	1 liter	BP1333-1
	20 liter	BP1333-20
Tris Borate EDTA, 1X Solution (TBE)	20 liter	BP2430-20
Tris Acetate EDTA, 50X Solution (TAE)	500 ml	BP1332-500
	1 liter	BP1332-1
	20 liter	BP1332-20
Tris Acetate EDTA, 1X Solution (TAE)	20 liter	BP2434-20

Accessories

Item	Catalog No.
Drip tray (for small incubator)	13-247-17
Drip tray (for medium/dual incubator)	13-247-27
Under cabinet mounting bracket (for small incubator)	13-247-18
Stacking plate (for small incubator)	13-247-19
Stacking plate (for medium incubator)	13-247-29
Rocker assembly	13-247-28
Hybridization tubes, 35 mm ID x 300 mm long (available in pack of 2 or case of 6)	13-247-300
Hybridization tubes, 35 mm ID x 150 mm long (available in pack of 2 or case of 6)	13-247-150
Hybridization tube cap O-ring	13-247-RNG
Rotisserie, 35 mm, 4/8 tubes	13-247-15
Rotisserie, 70 mm, 2 tubes	13-247-16
Rotisserie, 35 mm, 10/20 tubes	13-247-25
Rotisserie, 70 mm, 4 tubes	13-247-26

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